


Lymactin-NK

Anti HER2 monoclonal antibody

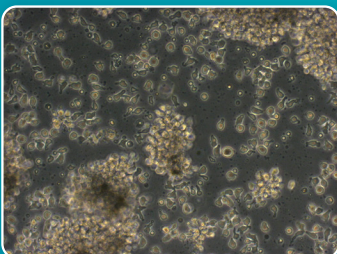

- Optimal NK cell activator for immune cell therapy research.
- Preservative free

Lymactin-NK

Lymactin-NK is anti HER2 monoclonal antibody. It can activate human and animal peripheral blood lymphocyte by immobilizing it on culture vessels, thus inducing NK proliferation.

- Used with ALyS505NK-AC, ALyS505NK-EX
- Preservative free ■ Conc.: 2.4mg/mL

Example of cell culture with Lymactin-NK

Fig.1 Cell Proliferation Curve


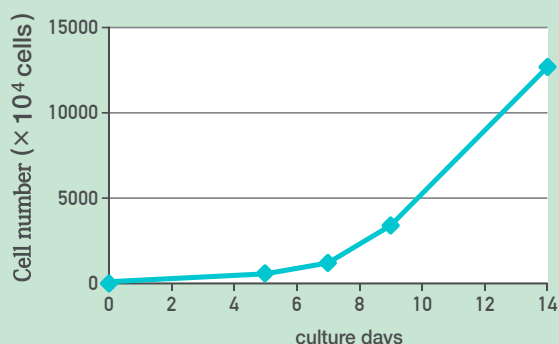


Fig.2 Change in Cell Surface Marker

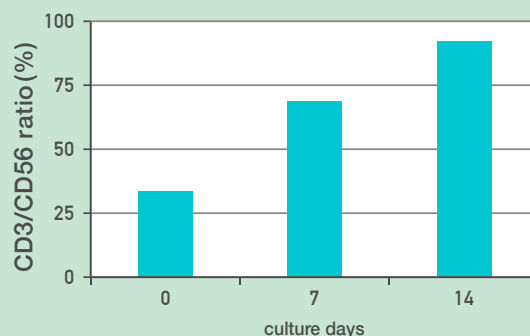

- Seeding density : 100×10⁴cells / mL • Culture scale : 12well plate → T225 Flask
- Medium : ALyS505NK AC1000 & EX1000 • Culture period : 14 days

Catalog# (CSTI / NIPRO)	Product Name	Description	Size (Packaging)
6002T01 / 87-025	Lymactin-NK	Anti HER2 monoclonal antibody (Preservative free)	2.4mg/mL (T)

(T) Tube

※This product is for research use only. It cannot be used for therapeutic or diagnostic purposes.



株式会社細胞科学研究所
宮城県仙台市青葉区西花苑1丁目16番16号
URL: <http://www.cstimedia.com>
E-mail: csti-info@ml.nipro.co.jp



ニプロ株式会社
大阪市北区本庄西3丁目9番3号

Official Distributor for North America
Iwai North America Inc.



Tel: 650-486-1541
orders@iwai-chem.com
www.iwaichem.com